



CELLIST_{TM} BASAL Medium (Model No. BASAL3P)

Name	Model number	Packaging	Contents	
CELLiST™ BASAL Medium	BASAL3P	1 L Aluminum Pouch	28.0 g	

Properties:

- Chemically-defined, animal origin-free, with no proteins or growth factors added.
- Does NOT contain hydrolysates, extracts or other undefined components.
- Does NOT contain proteins such as growth factors.
- Does NOT contain thymidine or hypoxanthine.
- Does NOT contain L-glutamine sources.
- Does NOT contain sodium bicarbonate.
- Contains poloxamer.
- Contains 10 g/L Glucose.

Storage conditions:

Before liquid preparation, store powder media in dark, refrigerated environment (2–8°C), away from high humidity. After liquid preparation, store in dark, refrigerated environment (2–8°C).

Instruction for preparation of liquid medium:

Table 1: Various parameters for preparation of 1 L of basal medium

	Powder weight	NaHCO₃ amount to be added	Weight of water to be added	pH*	Osmotic pressure*	Total solution weight	Specific gravity (Room temperature)
	28.0 g	1 9 a	986 mL	6.8-7.2	310-330	1010 g	1.01
28.0 g	1.8 g	(986 g)	0.0-7.2	mOsm/kg	1010 g	1.01	

^{*}Reference value

- 1. Prepare a suitable container and stir bar (magnetic bar).
- 2. Fill the container with about 90% volume (900 mL) of cell culture-grade water (room temperature).
- 3. The total amount of this pouch (28.0 g) should be added to the container. Place a small amount of cell culture-grade water in the pouch to wash the remaining product into the container.
- 4. Add 1.8 g of Sodium Bicarbonate.
- 5. Mix using magnetic stirrer for 20 minutes (until all powder is dissolved).
- 6. Add cell culture-grade water to final volume of 1 L and mix the media for 10 minutes. Volume adjustments can also be done by weighing (see table above).
- 7. Check pH to ensure a proper range of 6.8-7.2. If out of range, adjust pH using HCl or NaOH solutions.
- 8. Filter the medium in a biosafety cabinet, using a membrane filter with pore size of 0.2 to 0.22 μm in diameter.
- 9. Store in a refrigerated (2–8°C), dark environment until use.
- 10. Right before use, add L-glutamine or AminoStable™ to the solution (2-6 mM final concentration is recommended). If required, add growth factors such as insulin or IGF-I.

Use:

- This product is a cell culture medium used for research applications. Do not use it for any other purpose.
- For use in manufacturing, and for any other inquiries, please contact the following: